



---

## ImmEdge™ PEN

Catalog Number H-4000

**Description:** Hydrophobic barrier pen for immunohistochemistry and *in situ* hybridization. ImmEdge™ Pen provides a heat-stable water-repellent barrier that keeps reagents localized on tissue sections.

**Unit Size:** 2 pen set

**Storage Conditions:** Room temperature with pen tip down and tightly capped.

### INSTRUCTIONS FOR USE

1. Shake vigorously with lid on.
2. Depress tip to saturate, if necessary.
3. Circumscribe tissue section with the ImmEdge™ Pen.
4. Replace pen cap. Store pen tip down.
5. Allow ImmEdge™ Pen reagent to dry before immersing slides in aqueous solutions.

### DISPOSAL OF REAGENTS

Dispose in an approved landfill.

MSDS available online at: [www.vectorlabs.com](http://www.vectorlabs.com)

For Professional Use Only.

Authorized Representative / Exclusive Distributor to UK and Ireland:  
Vector Laboratories Ltd  
3 Accent Park  
Bakewell Road, Orton Southgate  
Peterborough, PE2 6XS, United Kingdom  
Tel: (01733) 237999 • Fax: (01733) 237119

A multi-lingual version of this data sheet can be found online at:-  
<http://www.vectorlabs.co.uk/3398/multi-language-instructions/>

Revised 03/16